

PRODUCT
INFORMATION

BRAND Reagent Reservoir 701460
and Liquid Handling Station

Reducing the residual volume

while working with the BRAND reagent reservoir 701460
in combination with a 50 µl multichannel liquid end

Abstract

Costly media, solutions, and samples are frequently used in everyday laboratory operations. In order to reduce costs, and minimize the environmental impact, it is of particular interest to work in the most resource-efficient way possible. The BRAND reagent reservoir 701460 is designed to minimize residual volume left behind during pipetting operations, especially for automated pipetting. As a result, it is ideally suited for automation. This technical note describes how the residual volume can be reduced even further while using the BRAND Liquid Handling Station.

Introduction

With the BRAND Liquid Handling Station (LHS), a wide variety of everyday laboratory pipetting tasks can be performed automatically, with accurate results. To prevent the occurrence of inaccuracies and errors, the LHS maintains a safety distance between the pipette tip and the bottom of the vessel, and between the pipette tip and the liquid surface. This ensures that no air is aspirated and no incorrect volumes are pipetted. Both safety distances result in a so-called “residual volume”, which remains in the vessel. Using the BRAND reagent reservoir with cover (Cat. No. 701460) and the 50 µl multichannel liquid end (50 µl MC LE) as an example, this technical note will demonstrate how the residual volume can be reduced through a simple change of the default settings in the LHS software.



50 µl multichannel
Liquid End

Material and method

To reduce the residual volume of the reagent reservoir with the LHS software, the reservoir is first added to the work table. By clicking on the labware, the options available for the labware appear on the right. The residual volume can be reduced in the “Details” tab after “use Custom Tolerances” has been activated (see Figure 1).

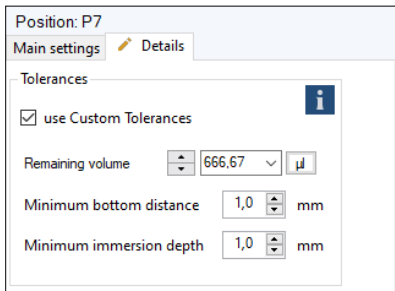


Figure 1
“Details” view of a labware

The safety distances mentioned previously are the so-called minimum bottom tolerance X_b and the minimum immersion depth X_i . X_b indicates the minimum distance between the pipette tip and the bottom of the vessel. The distance will not fall below this value. X_i indicates the minimum immersion depth of the pipette tip into the liquid. The sum of both safety distances is the liquid level that remains when the reservoir is emptied. This liquid level results in a residual volume V_r (Figure 2). X_b and X_i are each preset to 1 mm by default, resulting in a liquid level of 2 mm and, in the case of the reagent reservoir, a residual volume V_r of 666.67 µl with a total volume of 49 ml. To reduce V_r , the values X_b and X_i are alternately and successively reduced by 0.1 mm in the software. In the case of selected reservoir, this corresponds to a reduction of the

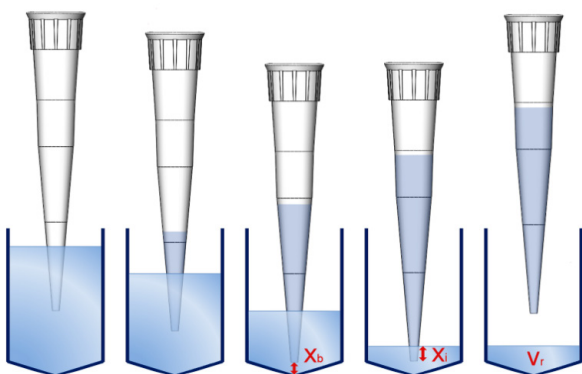


Figure 2
Graphical representation of the minimum bottom tolerance X_b , the minimum immersion depth X_i and the residual volume V_r .

residual volume V_r by 33.33 µl each time. Then, the 50 µl MC LE carries out a pipetting operation to remove the liquid up to the residual volume and to check whether a deterioration of the transfer can be detected. A deterioration manifests itself, for example, when air is aspirated or the pipette tip touches the bottom of the vessel.

In the transfer settings, the LHS is set to dispense above the reservoir so that the tips are more visible (Figure 3).

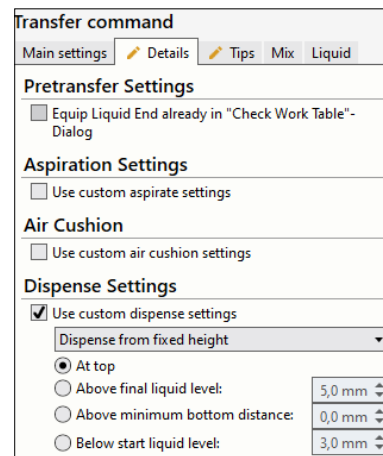


Figure 3
“Details” - view of a transfer.

In addition, prewetting of the pipette tips is deactivated (Figure 4), in order to prevent errors (see Results and discussion). Fully-deionized water (DI water) is used as the medium and the liquid type “Standard” is selected.

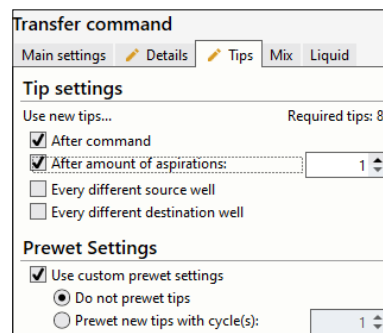


Figure 4
“Tips” - view of a transfer.

Results and discussion

As described above, the values for X_b and X_i were each reduced by 0.1 mm in succession, and as a result the residual volume V_r was reduced by 33.33 μl each time. Following this, a transfer removed the liquid up to the new residual volume and the pipetting quality was tested. The data is presented in the graph in Figure 5.

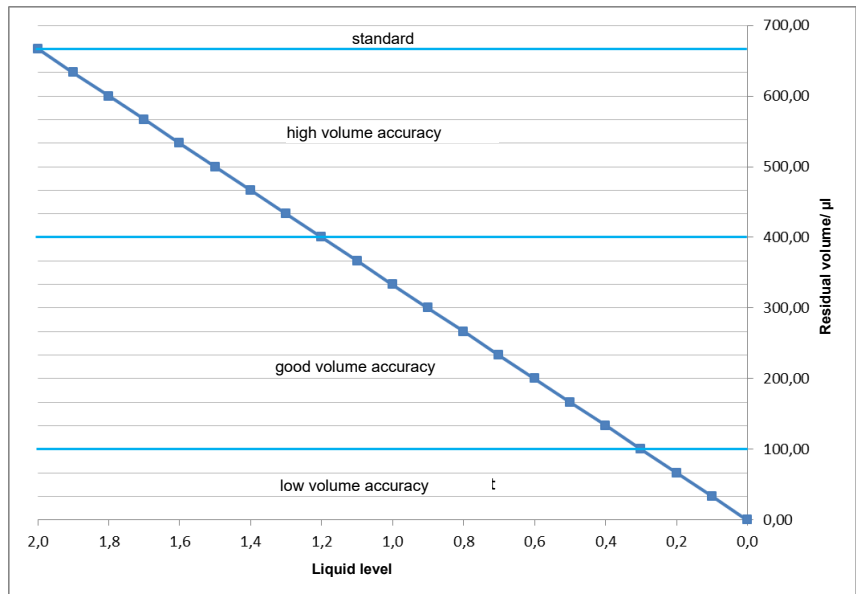


Figure 5
Graphical representation of the residual volume in μl plotted against the safety distance in mm. The different quality ranges can be seen.

Right at the beginning, it was apparent that the volume fill level on the left and right of the reservoir is higher than in the middle. This is probably due to adhesion forces with the vessel wall and depends on the liquid in use. It resulted in air being dispensed into the liquid because the tips were prewetted. The subsequent step of aspiration then drew in air along with the liquid. As a result, prewetting the tips was omitted for the entire experiment (Figure 4, right).

The pipetting operation was problem-free up to a residual volume V_r of 400.00 μl (liquid level = 1.2 mm) (see Figure 5). At a considerably lower residual volume, there was a risk of the tips touching the bottom of the reservoir. The potential contact between the tips and the bottom of the vessel is caused by the different tolerances of the liquid end, the

reservoir and by inserting the tips, and it illustrates the need for the safety distance X_b . As already mentioned, a liquid level of 1.2 mm in this case means that a minimum distance to the bottom X_b of 0.6 mm and a minimum immersion depth X_i of 0.6 mm were selected (compare with Figure 2, right). From a residual volume V_r of 66.67 μl or less (liquid level ≤ 0.2 mm), the volume was no longer pipetted correctly. Aspiration of air could be observed. This demonstrates that the residual volume V_r can be reduced to 100 μl (liquid level = 0.3 mm) without problems.

In general, however, when reducing the liquid level to between 0.3 and 1.1 mm (residual volume $V_r = 100.00 - 366.67 \mu\text{l}$), the dispensed volume should be checked.

Conclusion

Using the 4 x 40 ml reservoir 701460, the experiment demonstrated how the residual volume V_r is reduced by lowering the minimal immersion depth X_i and the minimum bottom tolerance X_b . As a result, the residual volume V_r for the 50 μl MC LE could be reduced from 666.67 μl to 400.00 μl or, as

the case may be, to 100.00 μl , depending on the accuracy required. This corresponds to a reduction of the default residual volume setting of approximately 40% (400.00 μl) or, as the case may be, approximately 85% (100.00 μl).



All information about our products can be found at shop.brand.de

BRAND GMBH + CO KG

P.O. Box 1155 | 97861 Wertheim | Germany

T +49 9342 808 0 | F +49 9342 808 98000 | info@brand.de | www.brand.de



BRAND. For lab. For life.®

BRAND®, BRAND. For lab. For life.®, as well as the BRAND figurative mark are registered trademarks or trademarks of BRAND GMBH + CO KG, Germany. All other trademarks mentioned or depicted here are the property of the respective owners.

Our technical literature is intended to inform and advise our customers. However, the validity of general empirical values, and of results obtained under test conditions, for specific applications depends on many factors beyond our control. Please appreciate, therefore, that no claims can be derived from our advice. The user is responsible for checking the appropriateness of the product for any particular application.

California Residents: For more information concerning California Proposition 65, please refer to www.brand.de/calprop65.

Subject to technical modification without notice. Errors excepted.

© 2021 BRAND GMBH + CO KG | Printed in Germany | 0621



Find accessories and replacement parts, user manuals, test instructions (SOP) and product videos at shop.brand.de



Further information on products and applications can be found on our YouTube channel: [mylabBRAND](https://www.youtube.com/mylabBRAND)

BRAND (Shanghai) Trading Co., Ltd.
Shanghai, China

Tel.: +86 21 6422 2318
info@brand.com.cn
www.brand.cn.com

BRAND Scientific Equipment Pvt. Ltd.
Mumbai, India

Tel.: +91 22 42957790
customersupport@brand.co.in
www.brand.co.in

BrandTech® Scientific, Inc.
Essex, CT. United States of America

Tel.: +1 860 767 2562
info@brandtech.com
www.brandtech.com